



Instructions for Use

- 1) Shake stick to end of bag.
- 2) Tear bag open.
- 3) Squeeze bag to open. Aseptically grasp the stick above the thumb-stop line to remove the sponge.
- 4) Aseptically swab across the entire sampling surface.
- 5) Turn sponge over. Change direction 90°. Aseptically swab the same sampling surface.
- 6) Aseptically place sponge into bag up to the thumb-stop.
- 7) Hold sponge in place inside bag. Bend stick to break. Allow sponge to drop in bag. Discard stick.
- 8) Fold bag to close.
- 9) Fold ends of blue wire inward.

Disposal: Follow established laboratory practices.



Storage Symbol: See chart.